**5 MEGA  
PIXELS**

**3.5"**

**LED  
ILLUMINATION**

**LI-ION  
BATTERY**

**500x  
ZOOOOM**

**USB 2.0**

**microSD**

# Levenhuk DTX 500 LCD

## Digital Microscope

User Guide

EN CZ PL RU

**levenhuk®**  
Zoom&Joy

# Levenhuk DTX 500 LCD

EN Digital microscope

Never, under any circumstances, look directly at the Sun, another bright source of light or at a laser through this device, as this may cause PERMANENT RETINAL DAMAGE and may lead to BLINDNESS.

## The kit includes

Microscope

Power cable

Cleaning cloth

Levenhuk DTX installation CD

User Guide


Rechargeable battery

USB cable


RCA cable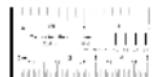

Calibration scale

# Microscope parts

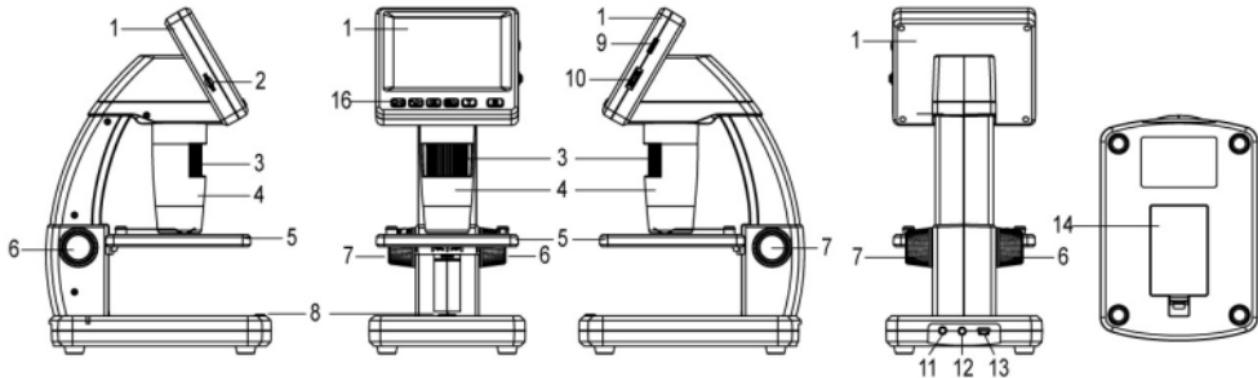- 1. LCD display
- 2. MicroSD card slot
- 3. Fine focus knob
- 4. Microscope optical tube
- 5. Stage
- 6. Coarse focus knob
- 7. Focus lock knob
- 8. Shutter button

- 9. Brightness adjustment wheel
- 10. Magnification adjustment wheel
- 11. Power cable socket
- 12. RCA socket
- 13. USB socket
- 14. Battery compartment
- 15. Battery status indicator
- 16. Function keys

# Microscope assembly

## Installing the rechargeable battery


Open the battery compartment at the bottom of the microscope base. Insert the battery into the battery compartment matching the indicated polarity.

*Note: the battery has to be fully charged before installation  
(refer to Charging the battery below).*

## Inserting the MicroSD card


A MicroSD card (purchased separately) is required to save captured images and recorded videos. Turn the microscope off and insert the MicroSD card into the corresponding slot on the microscope body until it locks in place with a click.

*Note: do not apply excessive force when inserting the MicroSD card.  
Try rotating the card, if it does not lock in place with ease. Format the card before using it.*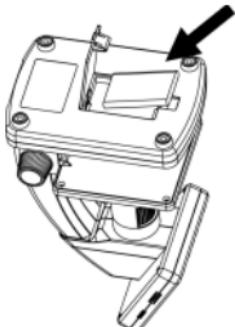

## Microscope display interface

-  Image capture mode icon
-  Sequence capture mode icon
- 00013/01618** Captured image counter/Total number of images that may be captured (based on the current resolution)
-  MicroSD status icon
- 5.0M** Current image resolution
-  Battery status indicator

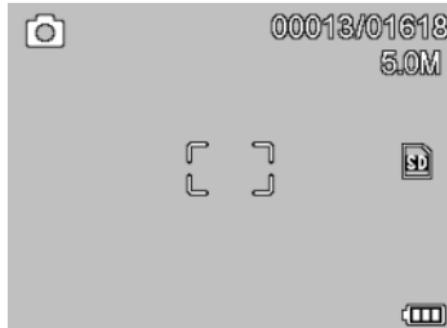*Note: interface icons may be hidden by pressing the magnification adjustment wheel (9).*

### Turning the microscope on

Press the power button (16).

### Turning the microscope off

Press the power button (16) again (refer to Power saving for more information).

# Using the microscope

## Capturing an image

Turn the microscope on and place a specimen on the stage. Adjust the illumination brightness, so that the specimen is evenly lit. Change magnification and focus your view, if necessary, then lock the results in place with the focus lock knob (7). Press the shutter button (8).

## Recording a video

Press the **◀** key (16) to switch to video mode. The icon in the upper left corner of the screen will change in confirmation. Press the shutter button (8) to start recording and press it again to stop.

## Playing recorded videos

Press the **◀** key (16) to switch to playback mode. The icon in the upper left corner of the screen will change in confirmation. Select the desired video clip with scroll keys (16) and press **OK** to play it.

## Deleting files

When browsing saved files, press the **Wrench** key (16), select **Delete...** press **OK** to confirm selection (you can choose to delete everything or just the selected file from the displayed menu).

## Protecting files

When browsing saved files, press the **Wrench** key (16), select **Protect** press **OK** confirm selection (you can choose to protect everything or just the selected file from the displayed menu).

## **Slideshow**

When browsing saved files, press the **Wrench** key (16), select **Slide Show** and press **OK** to confirm selection.

## **Interface color mode**

When browsing saved files, press the **▲** key to switch between existing interface color modes.

## **Magnification**

Scroll the magnification adjustment wheel (10) to zoom your view in or out.

*Note: magnification range of Levenhuk DTX 500 LCD is 1x-4x.*

## **Transferring files to a computer**

Connect the microscope to your computer with a USB cable and, when the AutoPlay window appears, choose to use the microscope as an external storage device.

## **Connecting the microscope to a computer**

Connect the microscope to your computer with a USB cable and, when the AutoPlay window appears, choose to use the microscope as a camera. Run the PortableCapture Pro application (refer to PortableCapture Pro software section for more information).

*Note: function keys are disabled when the microscope is connected to a computer as a camera. The microscope is operated through the PortableCapture Pro application. All the images and videos will be saved to the computer's hard drive.*

## **Charging the battery**

The battery has to be fully charged before turning the microscope on for the first time. Plug the power cable into the power cable socket on the microscope body (11) and connect it to a power source.

The battery status indicator (15) will turn red. Once the battery is fully charged, the indicator will go out. You can now disconnect the microscope from the power source and turn it on.

The fully-charged battery allows you to use the microscope for two hours. The illumination will grow dim and the image quality will decrease to inform you that the battery is low. Charge the battery, as described above.

## **Connecting the microscope to a TV-set**

Plug the RCA cable into the microscope and connect it to your TV-set (refer to the User Guide of your TV-set for additional information). Turn on the TV and select the correct channel from the corresponding menu of the TV-set (refer to the User Guide of your TV-set for additional information).

# Operating the microscope

## Main menu

Press the **Wrench** key (16) to open the main menu. From this menu you can set image resolution, adjust image quality and sharpness, as well as set parameters for sequence capture.

## Resolution

Select **Resolution** from the main menu choose the required image resolution. Press **OK** to confirm selection.

*Note:* model supports the following resolutions: VGA / 1.3M / 3M / 5M / 9M / 12M.

## Sequence capture

Select **Timer** from the main menu and set the total number of images to be captured in the **Frames** dialog window. Confirm your selection by pressing **OK**. A new dialog window will appear where you can set the interval between images using the **▲** and **▼** keys. Press **OK** to confirm and **◀** to return to capture mode. Press the shutter button to start capturing images. The icon in the upper left corner will change in confirmation.

Once the set number of images has been captured, the screen will return to standard capture mode. Press the shutter button at any time to stop sequence capture.

## **Quality**

Select **Quality** from the main menu and choose one of the existing settings. This will change the compression rate of captured images.

## **Sharpness**

Select **Sharpness** from the main menu and choose one of the existing settings.

## **Language**

Select **Language** from the main menu and choose one of the existing settings. Press **OK** to confirm selection.

## **Power saving**

Select **Power Save** from the main menu and set the idle time before the microscope automatically shuts down.

## **Display frequency**

Select **Frequency** from the main menu and choose one of the existing settings. Press **OK** to confirm selection.

## Date and time display

Select **Date Stamp** from the main menu and choose one of the existing settings. You can opt to display the date or both date and time. Press **OK** to confirm selection.

## TV connection

Select **TV Out** from the main menu and choose one of the existing settings. Press **OK** to confirm selection.

## Date and time

Select **Date/Time** from the main menu and set the current date and time. Press **OK** to confirm entry. You can also choose one of the following date formats: YY/MM/DD, DD/MM/YY or MM/DD/YY.

## Format SD card

Select **Format SD** from the main menu to format the installed memory card.

## Settings reset

Select **Sys. Reset** from the main menu to revert back to factory settings.

## Hardware version

Select **Version** from the main menu to check the current hardware version.

# PortableCapture Pro software

## Software installation

Insert the installation CD into your CD drive. The Main Menu window will be displayed. Select **PortableCapture software** to install the PortableCapture Pro application. Keep in mind that a system restart will be required during installation. The installation CD contains a copy of this User Guide, which you can read by clicking the corresponding entry in the Main Menu.

## First steps

Connect the microscope to your PC with a USB cable and run the PortableCapture Pro application. A pop-up message will warn you if the microscope is not connected to the PC.

## Toolbar icons

Capture an image

Start and stop  
video recording

Sequence capture (you can set  
the time to begin capturing  
images or videos, interval  
between them, video duration  
and the number of images  
or videos to be captured)

Full screen mode.  
To exit full screen mode,  
press **Escape** your keyboard  
or double-click anywhere  
on the screen

## Preview window toolbar

Save as

Copy to clipboard

Copy file

Previous image

Next image

Undo the last action

Repeat the last action

Draw

Text

Measurement

Calibration

## Capturing an image

Click the corresponding icon on the toolbar or select Photo from the Capture menu.

## Recording a video

Click the corresponding icon on the toolbar or select Video from the Capture menu.


## Sequence capture

Click the corresponding icon on the toolbar. A dialog window will appear where you can set a number of parameters for sequence capture.

Click the **Photo** radio button to set the time to begin capturing images, interval between them and the total number of images.

Click the **Video** radio button to set the time to begin recording videos, their duration, interval between them and the total number of clips.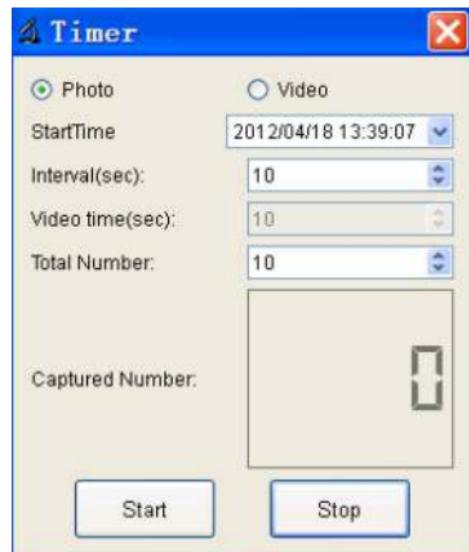



# Calibration and measurements

It is recommended that you calibrate the system before proceeding with measurements.

## Calibrating

Before calibrating the microscope, turn the grid display on. To do so, select **on** from **Options > XY Guide**. By default, the grid is visible. Place the calibration scale on the stage and focus your view. Make sure that the vertical axis of the grid is parallel to scale divisions, and capture an image.

Double-click the image thumbnail to open it in a separate window.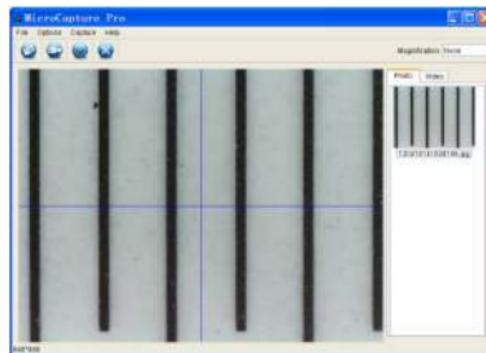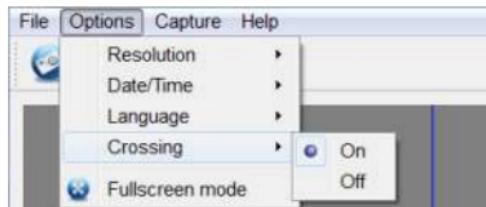


Click the corresponding icon on the toolbar to begin calibration. Using your mouse, choose two points on the image (you have to know the actual distance between these two points). Once you choose the second point, a dialog window will open, where you have to enter the known distance in the **Actual dimension** field.

The application will automatically calculate the magnification of the image. Click **OK** to close the dialog window. The calculated magnification will be displayed in the **Magnification** field.

**Note:** chosen points should form a horizontal line.

You can check calibration results by measuring the same distance with the **Any Angle Line** tool. Click the measurements icon, choose the **Any Angle Line** tool and draw a similar line on the image. If the measured distance is equal to the actual distance, calibration has been successful.

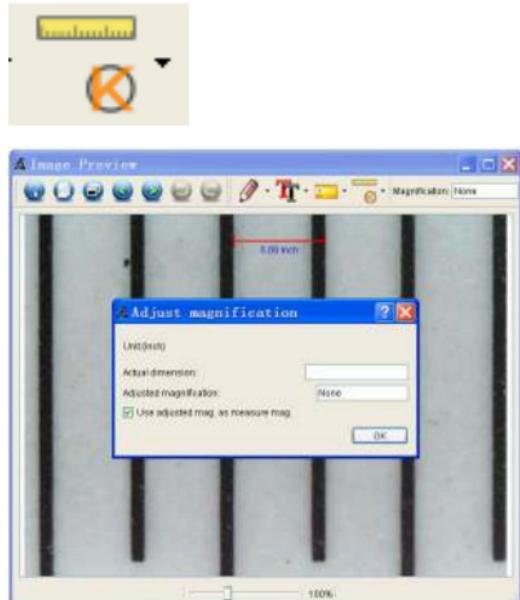This section explains all the measurement tools.

## Any Angle Line

Press and hold the left mouse button to start drawing a line.  
Release the left mouse button to complete the line.

## Continuous Line

Use this tool to draw a curve on the image and measure its length.

## Radius Circle

Draw a straight line on the image. The application will automatically calculate the radius, as well as circumference and area of the corresponding circle.

## Diameter Circle

Draw a straight line on the image. The application will automatically calculate the diameter, as well as circumference and area of the corresponding circle.

## Three Points Angle

Place three points on the image with your mouse. The application will automatically calculate the angle value.

# Recalibration

The calibration process has to be repeated if the magnification or focus has been changed during observations.

Click the corresponding icon on the toolbar again and select **Reset picture magnification** from **Reset magnification**.

Repeat the calibration process, as described above.

Change the magnification and capture a few images. Double-click one of the thumbnails to open the resulting image in a preview window. Click the calibration icon on the toolbar and select **Set picture magnification** from **Set magnification**. Enter the current magnification in the dialog window that appears and click **OK**.


If you want to use the same magnification when capturing images in the future, click the calibration icon on the toolbar and select **Set capture magnification** from **Set magnification**.


## Draw and text

PortableCapture Pro allows you to add drawings and textboxes to your images.

Open an image in a preview window and click the Pencil icon on the toolbar. Select one of the tools from the drop-down menu and draw something on the image.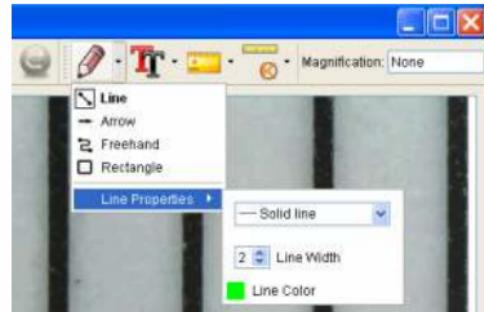

By clicking the TT icon on the toolbar, you can edit font type and color to be used in textboxes.

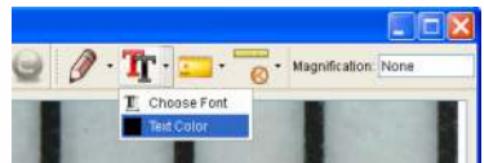# Working with the application for Mac OS users

Browse to the **mac** folder on the installation CD and copy **PortableCapture Pro** file to the desktop.

Operating the microscope  
Use a USB cable to connect the microscope to your computer.  
Run the PortableCapture Pro application by double-clicking its icon.

A magnified image of the observed specimen may be seen in the main window of the application.

From the **Options** menu, select **Preview Size** to set the image resolution. Keep in mind that the preview window cannot be resized (standard resolution is 640x480). The image resolution is shown in the bottom left corner of the preview window.

Click the corresponding icon on the toolbar or select **Photo** from the **Capture** menu to capture an image. An image thumbnail will appear on the right side of the main window.

Click the corresponding icon on the toolbar or select **Video** from the **Capture** menu to start recording a video. Editing images and videos  
Image and video thumbnails are displayed on the right side of the main window of the application.

*Note: during recording, the camcorder icon on the toolbar will turn red. Click the icon again to stop recording. A video thumbnail will appear on the right side of the main window.*

Double-click an image thumbnail to open it in a separate preview window. Editing images with PortableCapture Pro application is similar to working with other applications in Mac OS.

## Care and maintenance

Never, under any circumstances, look directly at the Sun, another bright source of light or at a laser through this device, as this may cause PERMANENT RETINAL DAMAGE and may lead to BLINDNESS. White LEDs are very bright; do not look at them as this may cause permanent damage to your vision. Take necessary precautions when using the device with children or others who have not read or who do not fully understand these instructions. After unpacking your microscope and before using it for the first time check for integrity and durability of every component and connection. Do not try to disassemble the device on your own for any reason. For repairs and cleaning of any kind, please contact your local specialized service center. Protect the device from sudden impact and excessive mechanical force. Do not apply excessive pressure when adjusting focus. Do not overtighten the locking screws. Do not touch the optical surfaces with your fingers. To clean the device exterior, use only special cleaning wipes and special optics cleaning tools from Levenhuk. Abrasive particles, such as sand, should not be wiped off lenses, but instead blown off or brushed away with a soft brush. Do not use the device for lengthy periods of time, or leave it unattended in direct sunlight. Keep the device away from water and high humidity. Be careful during your observations, always replace the dust cover after you are finished with observations to protect the device from dust and stains. If you are not using your microscope for extended periods of time, store the objective lenses and eyepieces separately from the microscope. Store the device in a dry, cool place away from hazardous acids and other chemicals, away from heaters, open fire and other sources of high temperatures. To avoid data loss or equipment damage, always close the application first, and then disconnect the microscope from your PC. Wait until the illumination goes off completely, and disconnect the microscope. Seek medical advice immediately if a small part or a battery is swallowed.

### Battery safety instructions

Always purchase the correct size and grade of battery most suitable for the intended use. Always replace the whole set of batteries at one time; taking care not to mix old and new ones, or batteries of different types. Clean the battery contacts and also those of the device prior to battery installation. Make sure the batteries are installed correctly with regard to polarity (+ and -). Remove batteries from equipment that is not to be used for an extended period of time. Remove used batteries promptly. Never attempt to recharge primary batteries as this may cause leakage, fire, or explosion. Never short-circuit batteries as this may lead to high temperatures, leakage, or explosion. Never heat batteries in order to revive them. Remember to switch off devices after use. Keep batteries out of the reach of children, to avoid risk of ingestion, suffocation, or poisoning.

# Specifications

Image sensor	5 Mpx (up to 12 Mpx by interpolation)
Magnification	20x-200x-500x
Available image resolutions	12 Mpx, 8 Mpx, 5 Mpx, 3 Mpx, 1.3 Mpx, VGA*
Photo/video format	.jpeg/*.avi
Memory card	microSD up to 32Gb (not included)
Objective lens	High-quality lens
Display	3.5" TFT-display, 4:3 ratio
Illumination	8-LED system with variable brightness
Power source	Li-ion battery (3.7 V/1050 mA·h) Operation time: 2 hours, charging time: 2 hours
Frame rate	30 fps
Power cable	100-240 V, 50/60 Hz; 5 V, 1 A output
OS language	English, German, French, Spanish, Italian, Portuguese, Japanese, Chinese
Software language	English, German, French, Spanish, Russian, Italian, Portuguese, Polish, Dutch, Japanese, Korean, Chinese
Dimensions (HxWxD)	227 x 104 x 149 mm
Weight	760 g (with the battery)

Levenhuk reserves the right to modify or discontinue any product without prior notice.

## System requirements

Operating system: Windows XP SP2/Vista/7/8, Mac 10.6-10.8,  
CPU: at least P4 1.8 GHz, RAM: 512 MB, interface: USB 2.0

# Levenhuk Limited Warranty

All Levenhuk telescopes, microscopes, binoculars and other optical products, except for accessories, carry a lifetime warranty against defects in materials and workmanship. Lifetime warranty is a guarantee on the lifetime of the product on the market. All Levenhuk accessories are warranted to be free of defects in materials and workmanship for six months from date of retail purchase. Levenhuk will repair or replace such product or part thereof which, upon inspection by Levenhuk, is found to be defective in materials or workmanship. As a condition to the obligation of Levenhuk to repair or replace such product, the product must be returned to Levenhuk together with proof of purchase satisfactory to Levenhuk.

This warranty does not cover consumable parts, such as bulbs (electrical, LED, halogen, energy-saving and other types of lamps), batteries (rechargeable and non-rechargeable), electrical consumables etc.

A Return Authorization (RA) Number must be obtained in advance of return. Contact the local Levenhuk branch to receive the RA number to be displayed on the outside of your shipping container. All returns must be accompanied by a written statement setting forth the name, address and telephone number of the owner, including a description of any claimed defects. Parts or products for which replacement is made will become the property of Levenhuk.

The customer will be responsible for all costs of transportation and insurance to and from Levenhuk or its authorized dealers and will be required to prepay such costs.

Levenhuk will use reasonable efforts to repair or replace any product covered by this warranty within thirty days of receipt. If a repair or replacement will require more than thirty days, Levenhuk will notify the customer accordingly. Levenhuk reserves the right to replace any product that has been discontinued from its product line with a new product of comparable value and function.

This warranty does not apply to any defects or damages resulting from alteration, modification, neglect, misuse, usage of improper power sources, damage in transportation, abuse, or any cause other than normal use, or to malfunction or deterioration due to normal wear. This warranty does not cover product that is lost, stolen, dropped, shattered, or any other type of damage or modifications to the original product caused by the customer. Levenhuk disclaims all warranties, express or implied, whether of merchantability or fitness for a particular use, except as expressly set forth herein. The sole obligation of Levenhuk under this limited warranty will be to repair or replace the covered product, in accordance with the terms set forth herein.

Levenhuk disclaims liability for any loss of profits, loss of information, or for any general, special, direct, indirect or consequential damages which may result from breach of any warranty, or arising out of the use or inability to use any Levenhuk product. Any warranties which are implied and which cannot be disclaimed will be limited in duration to a term of six months for accessories from the date of retail purchase.

Some states/provinces do not allow the exclusion or limitation of incidental or consequential damages, so the above limitations and exclusions may not apply to you. This warranty gives you specific legal rights, and you may have other rights which vary from state to state or province to province.

Levenhuk reserves the right to modify or discontinue any product without prior notice.

NOTE: This warranty is valid to USA and Canadian customers who have purchased this product from an authorized Levenhuk dealer in the USA or Canada. Warranty outside the USA or Canada is valid only to customers who purchased from an authorized Levenhuk dealer in the specific country or international distributor. Please contact them for any warranty service.

If warranty problems arise, or if you need assistance in using your product, contact the local Levenhuk branch:

Levenhuk Worldwide:  
USA: [www.levenhuk.com](http://www.levenhuk.com)  
Canada: [www.levenhuk.ca](http://www.levenhuk.ca)  
Czech Republic: [www.levenhuk.cz](http://www.levenhuk.cz)  
Germany: [de.levenhuk.com](http://de.levenhuk.com)  
Netherlands: [nl.levenhuk.com](http://nl.levenhuk.com)  
Poland: [pl.levenhuk.com](http://pl.levenhuk.com)  
Russia: [www.levenhuk.ru](http://www.levenhuk.ru)  
Ukraine: [www.levenhuk.ua](http://www.levenhuk.ua)  
EU: [www.levenhuk.eu](http://www.levenhuk.eu)

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The original Levenhuk cleaning accessories

# Levenhuk Cleaning Pen LP10

- Removes dust with a brush
- The soft tip is treated with a special cleaning fluid that removes greasy stains
- Does not damage optical coatings of the lenses
- Leaves no smudges or stains

[www.levenhuk.com](http://www.levenhuk.com)

Levenhuk, Inc. 1935 Brandon Court, Suite A-1, Chicago, IL 60139 USA  
Levenhuk® is a registered trademark of Levenhuk, Inc.  
© 2006-2015 Levenhuk, Inc. All rights reserved.  
20141111

**levenhuk**  
Zoom&Joy